

Product Information

PhasePrep™ BAC DNA Kit

Product Code NA 0100

Technical Bulletin Code MB-1035

TECHNICAL BULLETIN

Product Description

The PhasePrep™ BAC DNA Kit offers a scalable and cost-effective method for isolating large-molecular weight plasmids such as Bacterial Artificial Chromosomes (BAC) from recombinant *E. coli* cultures. The same kit can be used for preparations of four different sizes; sufficient reagents are provided for 300 micro, 180 mini, 30 midi, or 15 maxi preparations. Up to 2, 12, 50, or 100 µg of BAC DNA can be recovered from 5, 40, 250, or 500 ml of overnight recombinant *E. coli* culture, respectively. The purified BAC DNA contains very low levels of endotoxin (≤10 EU/µg DNA).

Recombinant *E. coli* culture is harvested by centrifugation and subjected to a modified alkaline-SDS lysis procedure. Nucleic acid is precipitated from the cleared lysate; residual RNA is removed by a short digestion at elevated temperature with an RNase cocktail. Endotoxins and other impurities are removed by simple temperature-mediated extraction and phase separation. Finally the BAC DNA is selectively precipitated from solution. The recovered BAC DNA is predominantly in its super-coiled form, free of RNA contamination, and ready for immediate use in downstream applications, such as sequencing, restriction digestion, cloning, and PCR[†].

Reagents Provided (Sufficient for 300 micro, 180 mini, 30 midi, or 15 maxi preparations)	Product Code	Quantity
Resuspension Solution	R 1149	2 x 200 ml
RNase A Solution (20 mg/ml)	R 6148	2 x 1.2 ml
Lysis Solution	L 1912	2 x 200 ml
Neutralization Solution	N 4659	2 x 200 ml
Elution Solution	E 5650	200 ml
RNase Cocktail	R 2151	250 µl
Endotoxin Removal Solution	E 4274	60 ml
Sodium Acetate Buffer Solution, 3M, pH 7.0	S 2404	25 ml
DNA Precipitation Solution	D 4564	200 ml

Storage

Store the kit at room temperature. Store the RNase Cocktail at -20 °C after opening the vial. The E4274 endotoxin removal solution may be stored at 2-8 °C for up to one month. For longer term, it is recommended to store the product at -20 °C."

Equipment and Reagents Required But Not Provided

- Ethanol, Product Code E 7148, E 7023, or 45,983-6
- Isopropanol, Product Code I 9516
- Centrifuge, capable of 15,000 x g, 2-8 °C
- Centrifuge tubes: 15 ml conical, Product Code C 3048; 50 ml Oak Ridge, Product Code T 2918
- Centrifuge bottles (250 ml & 500 ml), Product Codes B 1908 and B 2033
- Microcentrifuge
- Microcentrifuge tubes (1.5 ml & 2.0 ml), Product Codes T 9661 and T 2795
- 37 °C and 60 °C water baths

Precautions and Disclaimer

The PhasePrep BAC DNA Kit is for laboratory use only, not for drug, household, or other uses. Wear gloves, safety glasses, and suitable protective clothing when handling any reagent provided in the kit. See the Material Safety Data Sheet (MSDS).

Preparation Instructions

1. **Resuspension Solution:** Briefly centrifuge the tube containing the RNase A Solution. Add the RNase A Solution to the Resuspension Solution (one vial of RNase A Solution per bottle of Resuspension Solution), and mix thoroughly prior to initial use. Store the Resuspension Solution at 2-8 °C after adding the RNase A Solution.
2. **Lysis Solution:** If the SDS in the Lysis Solution precipitates during storage, heat the bottle at 37 °C until the SDS goes back into solution. Mix well and cool to room temperature before use.

3. **Neutralization Solution:** Chill on ice before use.
4. **Endotoxin Removal Solution:** Mix briefly and chill on ice before use.
5. **Heat water baths to 60 °C and 37 °C.**
6. **Centrifuge Speed:** To convert *g*-force to rpm or vice versa, refer to Appendix 1 for information.

Procedure

A. Micro Scale Preparation

1. **Harvest cells.** Pellet 3-5 ml of an overnight culture at 3,000-5,000 x *g* for 5-10 minutes. For best results, use a cell mass of 18 (cell mass = OD₆₀₀ x ml of culture; volume to use is equal to 18 divided by the absorbance of the culture at 600 nm). Remove the entire medium supernatant.

Note: The PhasePrep BAC DNA Kit is optimized for use with cultures grown in standard LB medium. For best results, inoculate the culture with a colony from a freshly streaked plate; grow the culture to OD₆₀₀ = 3-4. **Do not exceed a cell mass of 20.**

2. **Resuspend cells.** Add 250 µl of Resuspension Solution to the bacterial cell pellet. Make sure that the RNase A Solution has been added to the Resuspension Solution. Completely resuspend the bacterial pellet by pipetting up and down. Transfer the resuspended cells into a 1.5 ml microcentrifuge tube.
3. **Lyse cells.** Add 250 µl of Lysis Solution to the resuspended cells. Mix immediately by gently inverting the tube 4-6 times and incubate at room temperature for 5 minutes. **Do not vortex. Do not allow the lysis reaction to exceed 5 minutes.** Harsh mixing or prolonged lysis will shear the genomic DNA, resulting in genomic DNA contamination in the final preparation.
4. **Neutralize.** Add 250 µl of chilled Neutralization Solution to the lysate. Mix immediately and thoroughly by gently inverting the tube 6-8 times. **Do not vortex or shake**, as this will result in shearing the DNA. Incubate the tube on ice for 5 minutes. Centrifuge at maximum speed (~16,000 x *g*) for 5 minutes at 2-8 °C to pellet the cell debris. Carefully transfer the cleared lysate (supernatant) with a pipette tip to a clean microcentrifuge tube.

Note: To avoid taking floating particulates while transferring the supernatant, position the tip below the surface of the supernatant before aspirating.

5. **Precipitate nucleic acids.** Add 450 µl of isopropanol at room temperature to the cleared lysate. Mix thoroughly by gentle inversion and centrifuge at maximum speed for 20 minutes at 2-8 °C. Remove the supernatant and wash the pellet with 100 µl of 70% ethanol at room temperature. Centrifuge at maximum speed for 5 minutes at 2-8 °C, remove the supernatant by pipetting, and air dry the pellet briefly (5 minutes).
6. **Digest residual RNA.** Add 500 µl of Elution Solution to resuspend the pellet. Dilute the RNase Cocktail 10-fold by mixing 1 µl of the cocktail with 9 µl of Elution Solution. Add 1 µl of the diluted cocktail to the tube, mix briefly, and incubate in a 60 °C water bath for 5 minutes. Store the diluted RNase cocktail at 2-8 °C for later use (stable for at least three months).
7. **Adjust salt concentration.** Add 40 µl of Sodium Acetate Buffer Solution (3 M, pH 7.0) to the tube and mix briefly.

Note: To avoid shearing the BAC DNA, use large orifice pipette tips. Large orifice tips are also recommended for pipetting the viscous Endotoxin Removal Solution in the next step.

8. **Remove endotoxins and other impurities.**
 - a. Add 100 µl of Endotoxin Removal Solution and mix thoroughly by inversion. Chill the tube on ice for ≥ 5 minutes. Mix 1-2 times during the incubation. The solution will become light blue and clear.
 - b. Warm the tube in a 37 °C water bath for 5 minutes. The solution will become cloudy. Centrifuge at maximum speed for 3 minutes at room temperature to separate the phases. The clear upper phase contains the BAC DNA. The blue lower phase contains endotoxins and other impurities.
 - c. Carefully transfer the clear upper phase into a new 1.5 ml microcentrifuge tube. Discard the blue lower phase.
 - d. **Optional:** Repeat steps a, b, and c, then continue to step 9.
9. **Precipitate DNA.** Add 540 µl of room temperature DNA Precipitation Solution. Mix thoroughly by gentle inversion and centrifuge at maximum speed for 20 minutes at 4 °C. Remove

the supernatant and wash the pellet with 150 μ l of room temperature 70% ethanol. Centrifuge again for 5 minutes and pipette off the supernatant. Repeat the wash with 50 μ l of room temperature 70% ethanol. Remove all of the supernatant and air-dry the pellet briefly (5 minutes). Dissolve the DNA in 20-25 μ l of sterile deionized water (for sequencing) or Elution Solution. Rinse DNA from the side of the tube with gentle tapping. Use approximately one third of the preparation for a single sequencing reaction or restriction digestion.

Note: DNA aggregates may not be visible during the precipitation. After centrifugation the pellet may be loosely attached to the side of the centrifuge tube. Mark the tube for orientation relative to the centrifugal force prior to centrifugation. Carefully pipette off the supernatant from the side opposite the pellet to avoid the risk of aspirating the pellet.

B. Mini Scale Preparation

1. **Harvest cells.** Transfer 30-40 ml of an overnight culture to an Oak Ridge style centrifuge tube (capable of $\geq 15,000 \times g$) and centrifuge at 4,000-5,000 $\times g$ for 10 minutes. For best results, use a cell mass of 140 (see description of cell mass in the Micro Scale Preparation procedure). Remove all of the medium supernatant.

Note: The PhasePrep BAC DNA Kit is optimized for use with cultures grown in standard LB medium. For best results, inoculate the culture with a colony from a freshly streaked plate, and grow the culture to $OD_{600} = 3-4$. **Do not exceed a cell mass of 160.**

2. **Resuspend cells.** Add 2 ml of Resuspension Solution to the bacterial cell pellet. Make sure that the RNase A Solution has been added to the Resuspension Solution. Completely resuspend the bacterial pellet by pipetting up and down.
3. **Lyse cells.** Add 2 ml of Lysis Solution to the resuspended cells. Mix immediately by gently inverting the tube 4-6 times and incubate at room temperature for 5 minutes. **Do not vortex or shake. Do not allow the lysis reaction to exceed 5 minutes.** Harsh mixing or prolonged lysis will shear the genomic DNA, resulting in genomic DNA contamination in the final preparation.

4. **Neutralize.** Add 2 ml of chilled Neutralization Solution to the lysate. Mix immediately and thoroughly by gently inverting the tube 6-8 times. **Do not vortex or shake**, as this will result in shearing the DNA. Incubate the tube on ice for 5-10 minutes. Mix again during the ice incubation by gently inverting the tube 2-3 times. Centrifuge at 15,000 $\times g$ for 20 minutes at 2-8 $^{\circ}C$ to pellet the cell debris. Carefully transfer the cleared lysate (supernatant) with a pipette to a clean centrifuge tube.

Note: To avoid taking floating particulates while transferring the supernatant, position the pipette below the surface of the supernatant before aspirating.

5. **Precipitate nucleic acids.** Add 3.6 ml of room temperature isopropanol to the cleared lysate. Mix thoroughly by gentle inversion and centrifuge at 15,000 $\times g$ for 20 minutes at 2-8 $^{\circ}C$ (mark the tube for orientation relative to the centrifugal force before centrifugation). Pour off the supernatant and wash the pellet with 2 ml of room temperature 70% ethanol. Centrifuge at 15,000 $\times g$ for 5 minutes at 2-8 $^{\circ}C$ and pipette off the supernatant. Use a pipette tip to remove the last drop of liquid and air-dry the pellet briefly (5 minutes).
6. **Digest residual RNA.** Add 650 μ l of Elution Solution to resuspend the pellet. Rinse the pellet off the side of the tube with gentle swirling and dissolve the DNA for 5 minutes. Add 1 μ l of the RNase Cocktail to the tube, mix briefly, and incubate in a 60 $^{\circ}C$ water bath for 10 minutes. Rinse the side of the tube 1-2 times with gentle swirling during the digestion to dissolve any residual DNA pellet.
7. **Adjust salt concentration.** After digestion, centrifuge the tube briefly (1 minute) to collect the liquid. Transfer the solution to a 1.5 ml microcentrifuge tube. Add 50 μ l of Sodium Acetate Buffer Solution (3 M, pH 7.0) and mix briefly.

Note: The DNA solution may contain a small amount of insoluble particulates, which will be removed during the endotoxin removal steps. To avoid shearing the BAC DNA, use large orifice pipette tips. Large orifice tips are also recommended for pipetting the viscous Endotoxin Removal Solution in the next step.

8. **Remove endotoxins and other impurities.**
- Add 120 μ l of Endotoxin Removal Solution and mix thoroughly by inversion for 30 seconds. Chill the tube on ice for \geq 5 minutes. Mix 1-2 times during the incubation on ice. The solution will become light blue and clear.
 - Warm the tube in a 37 °C water bath for 5 minutes. The solution will become cloudy. Centrifuge at maximum speed (\sim 16,000 x g) for 3 minutes at room temperature to separate the phases. The clear upper phase contains the BAC DNA. The blue lower phase contains endotoxins and other impurities.
 - Carefully transfer the clear upper phase into a new 1.5 ml microcentrifuge tube. Discard the blue lower phase.
 - Repeat steps a and b, then continue to step 9.
9. **Precipitate DNA.** Transfer the clear upper phase into a 2 ml microcentrifuge tube. Add 700 μ l of room temperature DNA Precipitation Solution. Mix thoroughly by gentle inversion and centrifuge at maximum speed for 20 minutes at 4 °C. Remove the supernatant and wash the pellet with 500 μ l of room temperature 70% ethanol. Centrifuge again for 5 minutes and remove the supernatant. Repeat the wash with 100 μ l of 70% ethanol. Remove all of the supernatant and air-dry the pellet briefly (5 minutes). Dissolve the DNA in Elution Solution or sterile deionized water (for sequencing). Rinse the DNA off the side of the tube with gentle tapping.

Note: DNA aggregates may not be visible during the precipitation. After centrifugation, the pellet may be loosely attached to the side of the centrifuge tube. Mark the tube for orientation relative to the centrifugal force before centrifugation. Carefully pipette off the supernatant from the side opposite the pellet to avoid the risk of aspirating the pellet. Dissolve the DNA overnight in a refrigerator before quantitation.

C. Midi and Maxi Scale Preparations

- Harvest cells.** Transfer 150-250 ml (midi) or 300-500 ml (maxi) of an overnight culture to a centrifuge bottle; centrifuge at 4,000-5,000 x g for 10 minutes. For best results, use a cell mass of 750 for midi or 1500 for maxi (see description of cell mass in the Micro Scale Preparation procedure). Remove all of the medium supernatant.

Note: The PhasePrep BAC DNA Kit is optimized for use with cultures grown in standard LB medium. For best results, inoculate the culture with a colony from a freshly streaked plate, and grow the culture to $OD_{600} = 3-4$. **Do not exceed a cell mass of 1000 for a midi prep or 1750 for a maxi prep.**

- Resuspend cells.** Add 12 ml (midi) or 24 ml (maxi) of Resuspension Solution to the bacterial cell pellet. Make sure that the RNase A Solution has been added to the Resuspension Solution. Completely resuspend the bacterial pellet by shaking on a shaker or pipetting up and down.
- Lyse cells.** Add 12 ml (midi) or 24 ml (maxi) of Lysis Solution to the resuspended cells. Mix immediately by gently inverting the bottle 4-6 times and incubate at room temperature for 5 minutes. **Do not vortex or shake. Do not allow the lysis reaction to exceed 5 minutes.** Harsh mixing or prolonged lysis will shear the genomic DNA, resulting in genomic DNA contamination in the final preparation.
- Neutralize.** Add 12 ml (midi) or 24 ml (maxi) of chilled Neutralization Solution to the lysate. Mix immediately and thoroughly by gently inverting the bottle 8-10 times. **Do not vortex or shake**, as this will result in shearing the DNA. Incubate the bottle on ice for 10 minutes. Mix again during ice incubation by gently inverting the bottle 4-6 times. Transfer the lysate to a 250 ml centrifuge bottle (if the lysate is processed in a 500 ml bottle) and centrifuge at 15,000 x g for 30 minutes at 2-8 °C to pellet the cell debris. Carefully transfer the cleared lysate (supernatant) with a pipette to a 250-ml centrifuge bottle.

Note: To avoid taking floating particulates while transferring the supernatant, position the pipette below the surface of the supernatant before aspirating. If a large amount of cellular debris is carried over, re-centrifuge the supernatant for 10 minutes to remove the cellular debris. Alternatively, pass the cleared lysate through a fast flow filter paper prewetted with distilled water.

5. **Precipitate nucleic acids.** Add 22 ml (midi) or 44 ml (maxi) of room temperature isopropanol to the cleared lysate. Mix thoroughly by gentle inversion and centrifuge at 15,000 x *g* for 20 minutes at 4 °C (mark the bottle for orientation relative to the centrifugal force before centrifugation). Pour off the supernatant and wash the pellet with 10 ml of room temperature 70% ethanol. Centrifuge at 15,000 x *g* for 5 minutes at 4 °C and pipette off the supernatant. Use a pipette tip to remove the last drop of liquid and air-dry the pellet briefly (5 minutes).
 - c. Carefully transfer the clear upper phase into a new tube. Discard the blue lower phase.
 - d. Repeat steps a and b, then continue to step 9.

 6. **Digest residual RNA.** Add 5 ml (midi) or 10 ml (maxi) of Elution Solution to resuspend the pellet. Rinse the pellet off the side of the bottle with gentle swirling and dissolve the DNA for 5 minutes. Add 5 µl (midi) or 10 µl (maxi) of the RNase Cocktail to the bottle, mix briefly, and incubate in a 60 °C water bath for 15 minutes. Rinse the side of the bottle 1-2 times with gentle swirling during the digestion to dissolve any residual DNA pellet.

 7. **Adjust salt concentration.** After digestion, centrifuge the bottle briefly (1 minute) to collect the liquid. Transfer the solution to a 15 ml conical tube. Add 400 µl (midi) or 800 µl (maxi) of Sodium Acetate Buffer Solution (3 M, pH 7.0) and mix briefly.
- Note:** The DNA solution may contain a small amount of insoluble particulates, which will be removed during the endotoxin removal steps. To avoid shearing the BAC DNA, use large orifice pipette tips. Large orifice tips are also recommended for pipetting the viscous Endotoxin Removal Solution in the next step.
8. **Remove endotoxins and other impurities.**
 - a. Add 0.9 ml (midi) or 1.8 ml (maxi) of Endotoxin Removal Solution and mix thoroughly by inversion for 30 seconds. Chill the tube on ice for ≥ 10 minutes. Mix 1-2 times during the incubation on ice. The solution will become light blue and clear.
 - b. Warm the tube in a 37 °C water bath for 5 minutes. The solution will become cloudy. Centrifuge the tube at 3,000-4,000 x *g* for 5 minutes at room temperature to separate the phases (using a swinging bucket rotor). The clear upper phase contains the BAC DNA. The blue lower phase contains endotoxins and other impurities.

 9. **Precipitate DNA.** Transfer the clear upper phase into a 50-ml centrifuge tube. Add 5.4 ml (midi) or 10.8 ml (maxi) of room temperature DNA Precipitation Solution. Mix thoroughly by gentle inversion and centrifuge at 15,000 x *g* for 30 minutes at 4 °C. Remove the supernatant and wash the pellet with 5 ml (midi) or 10 ml (maxi) of room temperature 70% ethanol. Centrifuge again for 5 minutes and remove the supernatant. Repeat the wash with 2 ml of 70% ethanol. Remove all of the supernatant and air-dry the pellet for 5-10 minutes. Dissolve the DNA in Elution Solution or sterile deionized water (for sequencing). Rinse the DNA off the side of the tube with gentle swirling.
- Note:** DNA aggregates may not be visible during the precipitation. After centrifugation in a fixed angle rotor, the pellet may form a thin film along the centrifugal side of the tube and may be difficult to see. Mark the tube for orientation relative to the centrifugal force before centrifugation. Pipette off the supernatant from the opposite side of the pellet. Do not over-dry the pellet. Make sure to rinse any DNA off the side of the tube with gentle swirling while dissolving the DNA. Dissolve the DNA overnight in a refrigerator (tilt the tube to the side where the pellet is expected) before quantitation.

Results

Recovery and purity of BAC DNA preparations may be determined by spectrophotometric analysis. The ratio of absorbance at 260 nm and 280 nm (A_{260}/A_{280}) should be 1.7-1.9. The size and quality of DNA may be determined by restriction digestion and agarose gel electrophoresis or pulsed-field gel electrophoresis.

References

1. Birnboim, H. C., and Doly, J., A rapid alkaline extraction procedure for screening recombinant plasmid DNA. *Nucleic Acids Res.*, **7**, 1513-1522 (1979).
2. Birnboim, H. C., A rapid alkaline extraction method for the isolation of plasmid DNA. *Methods Enzymol.*, **100**, 243-255 (1983).
3. Cotton, M. *et al.*, Lipopolysaccharide is a frequent contaminant of plasmid DNA preparation and can be toxic to primary human cells in the presence of adenovirus. *Gene Therapy*, **1**, 239-246 (1994).

Troubleshooting Guide

Problem	Cause	Solution
Low DNA yield	Number of cells is insufficient	Inoculate the culture with a colony from a freshly streaked plate. For midi and maxi preps, grow a starter culture for 6-8 hr and inoculate an overnight culture with the starter culture in a 500 to 1000-fold dilution. Confirm that the cells were grown under optimal conditions (37 °C, 280-300 rpm) in a vessel at least four times larger than the volume of the medium. Grow the overnight cultures to OD ₆₀₀ = 3-4.
	Culture is too old	Streak a new plate from stock.
	Antibiotic activity is insufficient	Confirm that the medium contains the appropriate selective antibiotic. Use a sufficient amount of fresh antibiotic solution for growth of overnight cultures (for chloramphenicol, use 20 µg/ml for an overnight culture). Most antibiotic solutions are light sensitive and degrade during long term storage at 2-8 °C.
Low DNA yield (continued)	Lysis reaction is incomplete	Reduce the cell density or the culture volume, or increase volumes of Resuspension Solution, Lysis Solution, and Neutralization Solution proportionally. Keep the Lysis Solution bottle tightly capped between uses to prevent acidification. Check Lysis Solution for SDS precipitation; redissolve SDS by warming at 37 °C and allow to cool to room temperature before use.
	Lysate is not mixed appropriately with Neutralization Solution	After adding the prechilled Neutralization Solution, mix immediately but gently, by inversion until a fluffy egg-white material forms. Mix again during the incubation on ice.
	DNA pellet is lost or not rinsed off from the side of the tube	Mark the tube for orientation before centrifuging to help locate the pellet. Pipette off the supernatant from the opposite side of the pellet; avoid disrupting the pellet. Rinse the DNA off the side of the tube by gentle tapping or swirling when re-dissolving the DNA pellet.
Sheared BAC DNA	Alkaline lysis is prolonged	Do not allow the lysis reaction to exceed 5 minutes.
	Excessive mixing causes shearing	Mix gently by inversion during lysis and neutralization. Do not shake or vortex.
	Shearing occurs during reconstitution	Reconstitute the DNA gently. Avoid vigorous pipetting. Use large orifice pipette tips for handling the BAC DNA solution.
Poor performance in downstream applications	The final BAC DNA eluate contains too much salt or protein	Perform DNA precipitation with room temperature isopropanol in Step 5 and room temperature DNA Precipitation Solution in Step 9. Do not perform the precipitation at -20 °C. Wash the pellet twice in the final step with room temperature 70% ethanol and remove all of the liquid with pipette tips.
	EDTA is interfering by chelating cofactors	Precipitate the desired amount of BAC DNA and reconstitute in sterile deionized water for sequencing or other enzymatic reactions.
Absorbance of the final product does not match the actual quantity of BAC DNA	BAC DNA is contaminated with RNA	Confirm that the RNase A Solution was added to the Resuspension Solution prior to first time use. Store Resuspension Solution at 2-8 °C after adding the RNase A Solution. RNase A may degrade due to high temperatures (> 65 °C) or prolonged storage at room temperature. Keep the RNase Cocktail at -20 °C for long term storage.
	BAC DNA is contaminated with chromosomal DNA	Do not vortex or shake during lysis or neutralization. Do not allow the lysis reaction to exceed 5 minutes. Do not use cultures that have grown for over 24 hours or are in the cell death phase. Do not freeze and thaw bacterial cell pellets.
	DNA is not completely redissolved	Be sure that the DNA is completely redissolved before quantitation.

Appendix 1. Conversion of Centrifugal Force (in units of *g*) to RPM for Common Rotors

Centrifuge	Rotor	Type*	Radius (cm)	RPM at 3,000 x <i>g</i>	RPM at 5,000 x <i>g</i>	RPM at 15,000 x <i>g</i>
Beckman						
Allegra 6	GH-3.8	SB	20.4	3,631	4,688	N/A
Allegra 21(R)	S4180	SB	16.1	4,081	5,268	N/A
Allegra 64	F0485	FA	9.0	**N/A	N/A	12,211
	F0685	FA	9.7	N/A	N/A	11,764
TJ-25	TS-5.1-500	SB	19.0	3,756	4,849	N/A
	TA-10-250	FA	13.7	N/A	N/A	9,901
<i>Rotors for older Beckman centrifuges</i>	JA-10	FA	15.8	N/A	N/A	9,215
	JA-14	FA	13.7	N/A	N/A	9,896
	JA-20	FA	10.8	N/A	N/A	11,146
	JS-13	FA	14.0	N/A	N/A	9,790
IEC						
MP4(R)	215	SB	13.0	4,537	5,857	N/A
	224	SB	35.9	2,733	3,528	N/A
PR-7000M	966	SB	24.5	3,310	4,274	N/A
B22M	877	FA	12.6	N/A	N/A	10,318
Sorvall						
	HB-4	SB	14.7	4,277	5,522	N/A
	HB-6	SB	14.6	4,284	5,531	N/A
	HS-4	SB	17.2	3,948	5,097	N/A
	SH-80	SB	10.1	5,142	6,639	N/A
	GSA	FA	14.5	N/A	N/A	9,604
	SA-300	FA	9.7	N/A	N/A	11,784
	SA-600	FA	12.9	N/A	N/A	10,179
	SE-12	FA	9.3	N/A	N/A	11,997
	SL-50T	FA	10.7	N/A	N/A	11,203
	SS-34	FA	10.7	N/A	N/A	11,203
Eppendorf						
5410		12	5.8	N/A	N/A	15,652
5415C	F45-18-11	18	7.3	N/A	N/A	14,000
5415D&R	F45-24-11	24	8.3	N/A	N/A	13,155
5417C,D,&R	F45-30-11	30	9.5	N/A	N/A	12,279

*SB = swinging bucket; FA = fixed angle;

**N/A = not appropriate for application

See table for spin speeds in rpm for selected common centrifuges and rotors. The correct rpm for unlisted rotors can be calculated using the formula:

$$RPM = \sqrt{RCF / 1.118 \times 10^{-5} r}$$

where *RCF* = required gravitational acceleration (relative centrifugal force) in units of *g*; *r* = radius of the rotor in cm; *RPM* = the number of revolutions per minute required to achieve the necessary *g*-force.

† The PCR process is covered by patents owned by Hoffman-LaRoche, Inc.

Related Products	Product Code	Related Products	Product Code
Water, Molecular Biology Reagent	W 4502	Gel Loading Solution	G 2526
LB Broth, EZMix™	L 7658	λ DNA marker, <i>Hind</i> III, 125-23,130 bp	D 9780
LB Agar, EZMix™	L 7533	PulseMarker™ DNA marker, 0.1-200 kb	D 2291
Agarose for pulsed field electrophoresis	A 2929	Ethidium bromide, aqueous, 10 mg/ml	E 1510
TAE Buffer (10X)	T 9650	AutoPAGE™ Plus 4.5% Acrylamide Solution for Automated Sequencing	P 8977
TBE Buffer (10X)	T 4415		

FC 06/01

Sigma brand products are sold through Sigma-Aldrich, Inc.
 Sigma-Aldrich, Inc. warrants that its products conform to the information contained in this and other Sigma-Aldrich publications.
 Purchaser must determine the suitability of the product(s) for their particular use. Additional terms and conditions may apply.
 Please see reverse side of the invoice or packing slip.